

CONTERM SCIENTIFIC LTD

TECHNICAL MEMORANDUM

PRODUCT : GP SERIES 5

No : 0039

=====

=

FROM : Russell Kirkwood

DATE : 23/2/93

TO : ALL AGENTS

=====

=SUBJECT : S.E.L REPLACEMENT MOTORS (GP010).

A new type of motor is now fitted to the size 1, 2 & 3 GP SERIES 5 cabinets. This motor is supplied complete with a semi-integral rotor and adaptor mounting plate and has the advantage of being better matched to the required airflow characteristics of the smaller cabinet sizes.

This motor is **ONLY** suitable for the size 1, 2 & 3 cabinets and **MUST NOT** be fitted to the larger sizes as it will not generate sufficient airflow for correct operation.

This motor is an S.E.L shaded pole type and has our part number GP010.

Any new size 1,2 & 3 ex-factory cabinets will also be fitted with a specially baffled false back to enhance the operation of the S.E.L motor and will therefore no longer allow the older style full size motor (GP007) to be fitted.

If the new motor (GP010) is fitted as a replacement for the earlier motor (GP007) it will both fit and function **BUT** the airflow performance will **NOT** be as good as before.

All cabinets (Size 1, 2 & 3) supplied with an Appliance number 93xxx and later **SHOULD** be fitted with the new motor.

When ordering a replacement motor the customer should refer to the operating manual to determine which type is fitted to the cabinet.