

CONTHERM SCIENTIFIC LTD

TECHNICAL MEMORANDUM

PRODUCT : MOTORS

No : 0041

=====

=

FROM : Contherm Scientific Ltd

DATE : 10/9/93

TO : All Purchasers of replacement Motors

=====

=SUBJECT : Contherm Part Number: GP007 and P007.

PROBLEM:

Increase in motor length may prevent cabinet top cover from fitting in Contherm cabinets with top position circulating fans.

REMEDY:

Have a hole cut in lid 110mm diameter to allow fan motor body to protrude through.

We regret this remedy, however the increase in motor life will make the effort worthwhile.

Please contact Warren at our Help Desk on 0064 4 568-8034 for further assistance if required.

WARNING:

Do **NOT** remove aluminium spacer between motor and mount plate.