

CONTHERM *Scientific Limited*

TECHNICAL MEMORANDUM

PRODUCT :PRECISION

No : 0084

=====

=

FROM : Contherm Scientific Ltd

DATE : 1/7/2003

TO : ALL AGENTS

SUBJECT: REPLACING A RECTANGULAR GP FUSEHOLDER

The original rectangular GP Fuseholder is no longer available and must be replaced by a fuseholder adapter kit if it becomes necessary to replace the fuseholder.

The fuseholder kit consists of a tubular Schurter 230V 10A rated fuseholder and a black thermoplastic rectangular adapter to allow it to 'click in' the same hole as the original.

The wiring connects to the rear of the new fuseholder using the same quick connect spade terminal fittings as before.

A screwdriver is required to unscrew the fuse release to replace a fuse.

All future GP cabinets will be supplied with the new fuseholder arrangement.