

TECHNICAL MEMORANDUM

PRODUCT :GP

No : 0089a

---

FROM : Contherm Scientific Ltd

DATE : 29/6/2004 (Revised 31/01/2013)

---


TO : ALL AGENTS

---

**SUBJECT:** Replacing S.E.L motor with E.G.O Motor on GP cabinets

The S.E.L motor (GP1010/2010) is no longer available for the GP range of cabinets, it has been replaced by a new E.G.O motor (GP1000). The E.G.O motor and fan blade are very similar to the original. When replacing a S.E.L motor in the field a small kit is required, this will be automatically supplied whenever a GP1000/GP1010/2010 motor is ordered as a spare part. The kit consists of the following items:

1. GP1000 E.G.O motor complete with rotor which has been prewired with blue and brown motor wires of the same length (400mm) as the original motor. The motor adaptor ring has been modified to have a labelled earth stud fitted so that the original earth wire on the cabinet can be used with the replacement motor. Please ENSURE that the earth wire is connected to this terminal.
2. A copy of THIS techmemo.

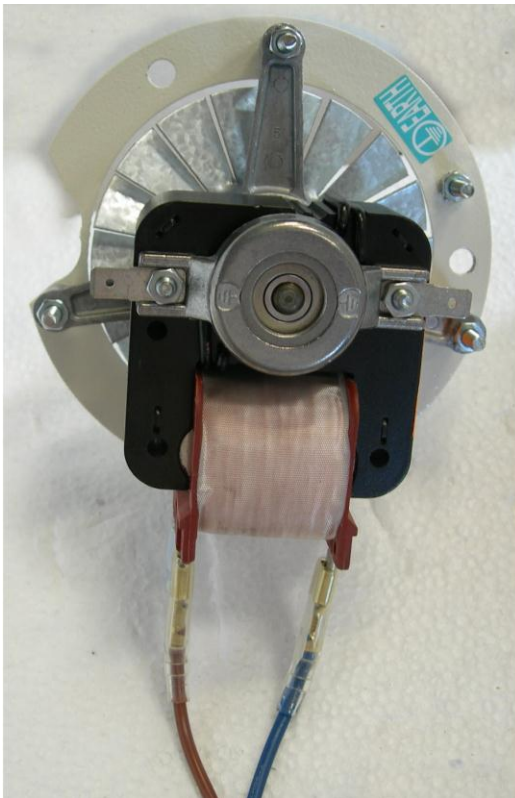

Connecting wires E.G.O motor (GP1000)



Fan blade